

DEXOLVE™ FOR IMPROVED PHARMACEUTICAL FORMULATIONS

STEP 1



Weigh in the following Dexolve™ amounts into 20 ml vials and prepare solutions with the given volume of distilled water:

Dexolve™*	Distilled water
3.0 g	7.0 mL
2.0 g	8.0 mL
1.0 g	9.0 mL
0.5 g	9.5 mL

*for accurate results take the water content of Dexolve™ into consideration

Use stirrer bar and magnetic stirrer.



DEXOLVE™ FOR IMPROVED PHARMACEUTICAL FORMULATIONS

STEP 2



- After the complete dissolution of Dexolve™, add ~50 mg or appropriate volume of your drug (candidate) to each vial. Should you be short of material, take smaller volume of the Dexolve™ solutions and dispense reduced amount of your substance, accordingly.
- Stir the resulting suspensions for 24 hours at room temperature. If your substance is sensitive, then cool your samples and protect them from light in the meantime.
- Observe the vials. If your substance completely dissolves upon stirring, dispense additional amount of your substance. Always ensure excess of material to be dissolved.



DEXOLVE™ FOR IMPROVED PHARMACEUTICAL FORMULATIONS

STEP 3



- When finished, filter the suspensions through PVDF syringe filters.
- Analyze the filtrate for your drug content.
- Establish relationship between the concentrations of Dexolve™ and the solubilized amounts of drug substance. Compare the data with the pure aqueous solubility of your substance.

In case you need technical help to facilitate the dissolution or to improve the solubilizing potency further, contact us!



Why use Dexolve™?

Possibilities

- Significant solubility enhancement (10 to 100,000 fold)
- Improvement of chemical stability
- Increased bioavailability, facilitated delivery
- Reduced aggregation
- Moderate irritation or reduced side-effects
- Maximized patient safety, complete renal elimination
- Enables formulation of water-insoluble APIs in all dosage forms
- Lower API doses can be achieved

APIs on the market or in development in formulations containing SBECD

- Voriconazole
- Carfilzomib
- Amiodarone
- Ziprasidone
- Maropitant (veterinary use)
- Posaconazole
- Carbamazepine
- Melphalan
- Topiramate
- Aripiprazole
- Clopidogrel
- Docetaxel

Several other nitrogen containing API bases are in various clinical phases.

DEXOLVE™ FOR IMPROVED PHARMACEUTICAL FORMULATIONS

COMPANY CONTACTS

CYCLOLAB CYCLODEXTRIN RESEARCH & DEVELOPMENT LABORATORY LTD.

Budapest, P.O. Box 435, H-1525 Hungary

Location: Illatos út 7., Budapest, H-1097- Hungary

Tel: (+36) 1-347-60-70

E-mail: dexolve@cyclolab.hu

Web: <http://www.cyclolab.hu>

CONTACT PERSON

Zoltán Kovács

Senior Business Development Specialist

kovacs@cyclolab.hu

(+36) 30-163-71-77

ASK FOR
A FREE
SAMPLE

